

Sartorius Stedim Biotech GmbH, 37070 Göttingen

To whom it may concern

July 12, 2023

Advisory Letter - Supplementary information for Minisart® HY | SRP Syringe Filters

Dear Valued Customer,

With this letter we would like to inform you of a possibly unclear description in the manual of Minisart® NML | HY | SRP Medical Device Syringe Filters for the following products:

Material number	Description
16596-----HYK	Minisart HY, hydrophobic PTFE, 0.2 µm, 26mm, sterile, 50 pack
16596-----HYQ	Minisart HY, hydrophobic PTFE, 0.2 µm, 26mm, non-sterile, 50 pack
16599-----HYQ	Minisart HY, hydrophobic PTFE, 0.2 µm, 26mm, non-sterile, 500 pack
16596-----ZSK	Minisart HY, hydrophobic PTFE, 0.2 µm, 26mm, sterile, 50 pack
17575-----ACK	Minisart SRP, hydrophobic PTFE, 0.2 µm, 25mm, sterile, 50 pack
17575-----ZSK	Minisart SRP, hydrophobic PTFE, 0.2 µm, 25mm, sterile, 50 pack

The current description in the manual indicates several possible applications. Please be informed that sterile filtration capacities of oily solutions and liquids are depending on the solution properties. This applies to all Minisart® HY | SRP syringe filters listed above, regardless of the manufacturing date and batch. **Please be advised that the sterile filtration capacities in your specific application for oily solutions and liquids must be tested and validated by the user according to the needs of the specific application.**

We thank you for your trust and continued use of Sartorius products. If you have any questions, please contact labfiltrationpm@sartorius.com.

Kind Regards

Adam Green
Head of Product Management
LPS Reagents & Consumables

Dr. Anna Vreemann
Site Quality Manager
Operations SSB Göttingen

Sartorius Stedim Biotech GmbH
August-Spindler-Straße 11
37079 Göttingen, Germany
Phone +49 551 308 0
Fax +49 551 308 3289
www.sartorius.com

Registered Office:
Göttingen
Local Court of Registration:
Amtsgericht Göttingen
HRB No. 200266

Managing Directors:
Michaela Pischke,
Dr. Helge Henning

Chairman of the Supervisory
Board:
Dr. René Fáber

Sartorius hereby expressly declares that this letter does not and shall not be deemed to constitute any acknowledgement, concession or admission that Sartorius and/or its Affiliates or suppliers have breached or violated any law, rule, contract or regulation, or an acknowledgement that you or your affiliates, customers, or any other third party may have any claims against Sartorius and/or its Affiliates relative to this subject matter or any that could have been asserted pursuant to these details.